

# Zyla sCMOS

## Introducing Zyla 4.2 PLUS

### 82% QE and Industry Fastest Speed




#### NEW ZYLA 4.2 PLUS

- 4.2 megapixel sCMOS
- 82% QE, optimized for all fluorophores
- 0.9 e<sup>-</sup> read noise
- 100 fps (53 fps USB 3.0)
- 33,000:1 dynamic range

#### ZYLA 5.5

- 5.5 megapixel sCMOS
- Rolling & True Global Shutter
- 1.2 e<sup>-</sup> read noise
- 100 fps (40 fps USB 3.0)
- 25,000:1 dynamic range

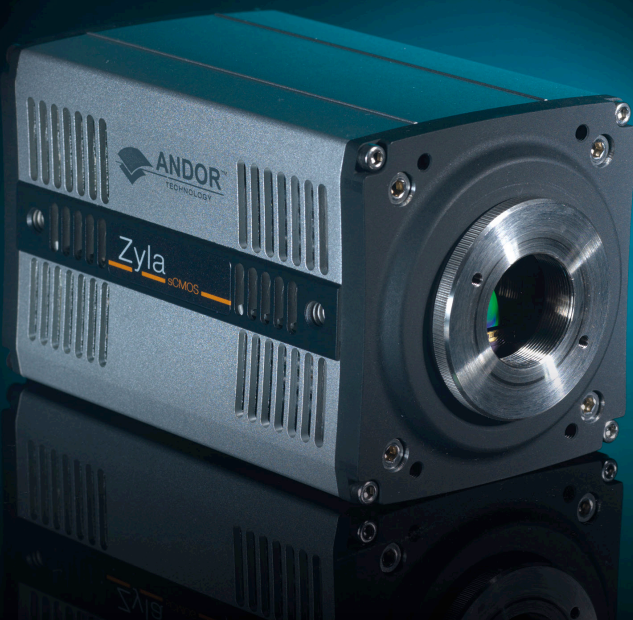


- ✓ QE Boosted to 82%
- ✓ Industry fastest USB 3.0 speeds
- ✓ LightScan PLUS
- ✓ >99.8 % Quantitative Linearity

# KEY INNOVATIONS OF THE NEW ZYLA 4.2 PLUS

## 1 QE Boosted to 82%

The latest-generation sCMOS sensor delivers a further 10% boost in QE, providing excellent broad coverage of the VIS/NIR region.

- ✓ Optimized for a broad range of fluorophores
- ✓ Reduced exposure times & faster frame rates
- ✓ Reduced phototoxicity / photobleaching
- ✓ Lower fluorophore concentrations – more accurate physiology

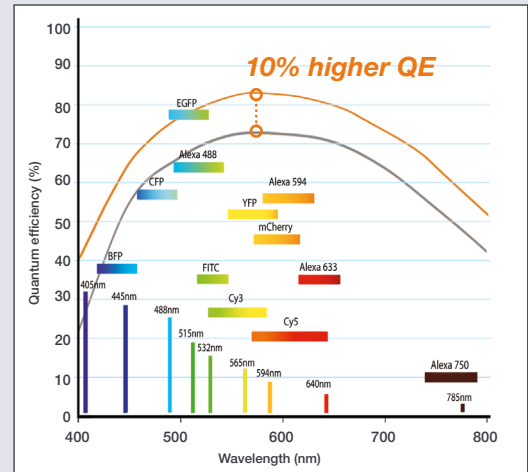

[Click here](#) to find out more about the Zyla 4.2 PLUS.



## 4 Application Modes

- LightScan PLUS** – Adapts the Rolling Shutter scan mode to applications such as Scanning LightSheet Microscopy and Line Scan Confocal
- FCS Mode** – Achieve up to 26,041 fps, ideal for Fluorescence Correlation Spectroscopy

## 2 Market Leading Speed

Superior data transfer efficiency and Zyla's unique 12-bit high speed mode combine to deliver an incredible 53 fps through super-convenient USB 3.0, up to 77% faster than competitor sCMOS cameras.

With this capability, speed is on tap to allow you to follow faster dynamic processes with improved temporal resolution. Opt for the Camera Link version to access up to a blistering 100 fps (full resolution).


- ✓ Industry fastest USB 3.0 frame rates
- ✓ Unique 12-bit mode for even faster speeds
- ✓ Up to 77% faster (full resolution) than competing sCMOS
- ✓ Unique ROI processing capability – Faster ROI speeds
- ✓ Follow faster temporal processes with a super-convenient USB 3.0 camera

Mode	Competing sCMOS USB 3.0	Zyla 4.2 PLUS USB 3.0
12-bit/ Full resolution	Not available	<b>53 fps</b>
16-bit/ Full resolution	30 fps	<b>40 fps</b>
12-bit/ 1024 x 1024	Not available	<b>200 fps</b>
16-bit/ 1024 x 1024	60 fps	<b>160 fps</b>

## 3 Market Leading Quantitative Linearity

Zyla 4.2 PLUS uses enhanced on-head intelligence to deliver market-leading linearity of > 99.8%, for unparalleled quantitative accuracy of measurement across the full dynamic range.

- ✓ Better than 99.8% linearity
- ✓ Increased quantitative accuracy

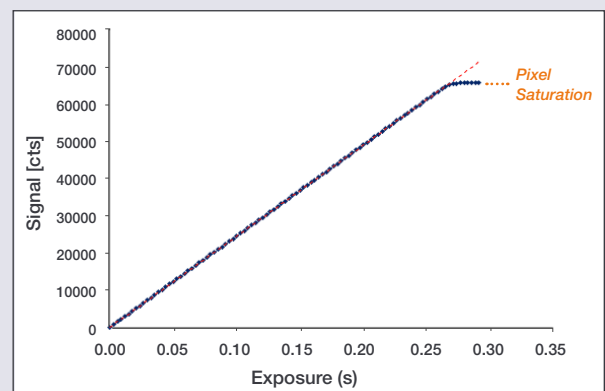

# FEATURES & BENEFITS




Andor's Zyla sCMOS camera platform offers high speed, high sensitivity and high resolution imaging performance. The remarkably light and compact, thermoelectrically-cooled design, integrates perfectly into both laboratory and OEM applications alike. Zyla is ideally suited to many cutting edge experiments that push the boundaries of speed and sensitivity.

The unprecedented value and flexibility of the Zyla means it is also re-defining the concept of a 'workhorse' camera, rapidly displacing interline CCDs as the gold standard microscope detector.

Feature	Benefit
<b>NEW</b> QE <sub>max</sub> boosted to 82%	Highest available photon capture efficiency across visible/NIR, optimized for all common fluorophores. Shorter exposures, reduced phototoxicity, lower dye concentrations for more accurate physiology.
~ 1 e <sup>-</sup> Read Noise	Noise floor down to 0.9e <sup>-</sup> . Lower detection limit than any CCD
<b>NEW</b> Market leading USB 3.0 speed	Superb USB 3.0 data transfer efficiency and Zyla's unique 12-bit high speed mode deliver up to 53 fps full resolution, 77% faster than competing sCMOS. Follow dynamic processes with improved temporal resolution.
100 fps (Camera Link)	Zyla offers '10-tap' Camera Link for maximum sustained frame rates. (Burst to 4GB on-head memory on Neo).
5.5 & 4.2 megapixel sensor formats and 6.5 μm pixels	Extremely sharp resolution over a 22 mm (Zyla 5.5) and 19 mm (Zyla 4.2 PLUS) diagonal field of view. Ideal for cell microscopy, astronomy and area scanning applications.
Rolling and Global shutter (Zyla 5.5)	Maximum exposure and readout flexibility across all applications. Global Shutter for 'interline CCD mode' freeze frame capture of fast moving/changing events.
Extended Dynamic Range	Unique 'dual gain amplifier' sensor architecture offering dynamic range of up to 33,000:1.
12-bit and 16-bit modes	12-bit Mode for smaller file size and absolute fastest frame rates through USB 3.0; 16-bit mode for full dynamic range
<b>NEW</b> Better than 99.8% linearity	Unparalleled quantitative measurement accuracy across the full dynamic range.
Very Low Fan Vibration	Implemented on both models. Designed with vibration sensitive experiments in mind, such as super-resolution microscopy
<b>NEW</b> LightScan PLUS with FlexiScan and CycleMax (Zyla 4.2)	Reduce background and improve contrast and resolution in scattering samples. Designed to allow users to maximise fluorescence signal and confocality concurrently in applications such as Scanned Light Sheet Microscopy and Line Scanning Confocal Microscopy
<b>NEW</b> FCS Mode	Fluorescence Correlation Spectroscopy requires the fastest possible speed from a minimal height ROI. Zyla 4.2 PLUS outputs a sustained 26,041 fps from a 2048(h) x 8(v) ROI.
Dark Noise Suppression (DNS) technology	Extremely competitive low darkcurrent of 0.14 e/pix/sec with fan cooling. Maintains low noise advantage across range of exposure conditions.
TE cooling to 0° C in up to 30 °C ambient	Ideal for OEM integration into enclosed systems.
Compact and Light	Ideal for integration into space restrictive set-ups. Ideal for OEM.
<b>NEW</b> GPU Express	Simplify and optimize data transfers from camera to Graphical Processing Unit (GPU) card to facilitate accelerated GPU processing as part of the acquisition pipeline.
Dynamic Baseline Clamp	Ensure quantitative stability
Hardware Timestamp	FPGA generated timestamp with 25ns accuracy.

## Zyla 5.5

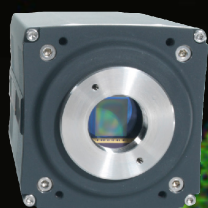

**Zyla 5.5** is truly unique in offering **both Rolling and true Global shutter** capability in one sensor. Global shutter offers 'snapshot' imaging capability, whereby all pixels in the area are exposed simultaneously, and is directly analogous to that which is available in interline CCDs. True Global shutter is only available through the '5T' (5-transistor) sensor design exploited in the Zyla 5.5 offering greater application flexibility and is ideal for tight synchronization with microscope peripheral devices such as z-stage or switchable light source.

Please see **page 6** for a further comparison of Rolling shutter and Global shutter modes.



# Zyla

## THE BIOLOGIST'S CHOICE

Zyla sCMOS has proven a superb camera choice for the biologist and microscopist. Many simply see the Zyla as an **amazing value**, superb price/performance 'workhorse' camera with which to replace their existing interline CCD and **upgrade** the performance of their fluorescence microscope. Others are driven by distinct **application performance criteria** that only sCMOS can answer.

### Quality, Throughput, Performance, Accessibility...

- **High Sensitivity & Wide Dynamic Range** – quantify very weak and very bright structures with one image.
- **Superb Image Quality** – high resolution and uniform backgrounds for publication quality imaging.
- **Capture Everything** – the larger field of view matches that of modern microscopes. Achieve better statistics and higher throughput in high content experiments.
- **Blazingly Fast** – more and more studies of cell processes require greater temporal resolution.
- **GPU Express** – for real time processing.
- **Ease of use** – designed to get you up and imaging with minimal fuss.
- **Flexible** – fast or slow, big or small, weak or bright... Zyla is adaptable for all of your imaging challenges.

### Example Areas of Application

#### Physiology / Ion Imaging

The fast frame rate and excellent sensitivity of Zyla is ideally suited to the particular needs of ion signalling microscopy. Zyla 4.2 PLUS offers superlative sensitivity at speed, but electrophysiology may require the Global Shutter exposure mode of Zyla 5.5 to ensure temporal correlation across the whole image.

#### Super Resolution Microscopy

The low vibration, high QE, low noise and speed capability of Zyla 4.2 PLUS (USB 3.0 and Camera Link) is well suited to the particular detection criteria of single molecule based 'STORM / PALM' approaches, and is used by some as an alternative to EMCCDs for this purpose. Note, this should be considered distinct from the general needs of single molecule microscopy, which are best served by back-illuminated EMCCD cameras (see Andor iXon EMCCD range). Capability to switch off interpolative filtering and provision of custom blemish maps. **GPU Express** for real time data processing.

#### High Content Screening

Zyla sCMOS yields markedly improved throughput and statistical validity of data in high content analysis. For example, a larger field of view results in analysis of more cells per image, wider dynamic range means a field of variable intensity cells can be quantified in only one acquisition, and higher sensitivity results in reduced acquisition times. **GPU Express** for real time data processing. For further information, view this article: [highcontentreview.com/scmos/](http://highcontentreview.com/scmos/)

#### Light Sheet Microscopy

Andor sCMOS cameras have been at the forefront of innovative Light Sheet Microscopy development and significantly the Zyla 4.2 PLUS is now equipped with **LightScan PLUS**.

**LightScan PLUS**, ensures the user has additional control and flexibility over the functionality of the rolling shutter scan mode. LightScan PLUS allows the user to scan their scanning light source from the top to the bottom of the sensor, or vice versa, in one continuous sweep. In addition to this, **FlexiScan** permits independent adjustment of the line readout speed, the exposure time and the pixel row height (slit width), allowing signal strength and confocality to be optimized *concurrently*. **CycleMax**, ensures the fastest frame rates can be achieved with no dead time and no need to reset the laser for each alternate frame. **GPU Express** for real time data processing.

#### Cell Motility

The motile cell is captured extremely well by the speed and resolution of the Zyla. Generally, the rolling shutter of Zyla 4.2 PLUS is suited, but care must be taken of distortive effects if the cell is moving particularly fast. For example, it has been noted that the Zyla 5.5 in global shutter mode was required to image motile sperm cells.



#### TIRF Microscopy

The Zyla's fine pixel resolution, great sensitivity, large field of view and fast imaging speed offers a superb choice of platform for following/tracking fast processes at the cell membrane. Multi-wavelength TIRF may benefit from Zyla 5.5 in global shutter.

### Upgrade your microscope performance using Zyla sCMOS

Zyla remains within the same price bracket as interline CCDs, yet offers remarkable performance improvements:

- 4x more pixels
- 5x more sensitive
- 10x more dynamic range
- 16x faster

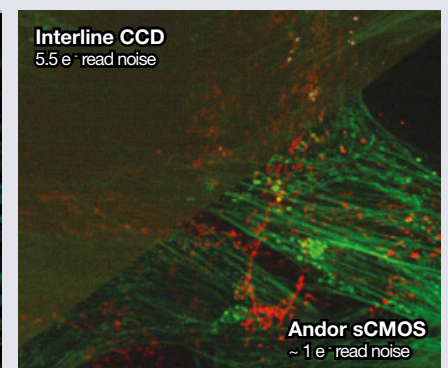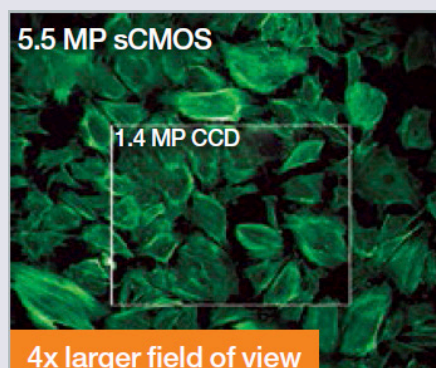

# Zyla

## THE PHYSICIST'S CHOICE

Zyla sCMOS has become a well established detector amongst physicists, biophysicists and astronomers, the advanced combination of speed, sensitivity and dynamic range enabling new ground to be broken.

### Performance & Adaptability

- **Dual Amplifier** – novel pixel architecture means you don't need to pre-select gain. Access lowest read noise and full well depth simultaneously.
- **1000 fps** – Access extremely fast frame rates through user definable Region of Interest control, suited to many applications within the physical sciences.
- **GPU Express** – for real time processing.
- **Global Shutter** – Zyla 5.5 offers this important mode that completely avoids spatial distortion, and ensures temporal correlation across all regions of the sensor. Achieve sub-microsecond inter-frame gaps in PIV applications.
- **Low darkcurrent** – low read noise is complimented by extremely competitive darkcurrent, also ensuring minimized hot pixel blemishes.
- **Cooling options** – standard Zyla 5.5 camera air cools to 0°C at up to +30°C ambient. Water cooled option available on request.
- **Blemish correction maps and advanced control** – Andor provide the capability to turn off/on blemish correction for those who prefer to perform this themselves. Bespoke blemish maps can also be provided.
- **Compact and Light** – the extremely small volume footprint of Zyla renders it adaptable to intricate optical set-ups.

### Example Areas of Application

#### Particle Imaging Velocimetry (PIV)

The true Global Shutter mode of Zyla 5.5 facilitates an inter-frame gap of down to 100 ns.

#### Lucky / Speckle Imaging

Zyla's fast frame rate and large field of view are ideal for this resolution enhancing technique. **GPU Express** for real time data processing.

#### Solar Astronomy

Fast frame rates, wide dynamic range and great linearity present a very formidable solution to the specific detector needs of next generation large solar telescopes.

#### Bose Einstein Condensation

The QE profile of Zyla is very good in the red/NIR region, ideal for BEC of Rb.

#### Adaptive Optics

Accessing > 1000 fps using ROIs renders the Zyla an ideal Wavefront detector. Use with data splitter to enable direct data access.

#### Fluorescence Correlation Spectroscopy

Superb temporal resolution from small ROIs are excellent for accurately measuring diffusion coefficients.

#### X-ray / Neutron Tomography

The Zyla can be readily lens coupled to scintillators and phosphors, presenting a high resolution, sensitive and fast solution for tomography.\*



sCMOS image of M81 and M82 in Ursa Major. Courtesy of Jin Ma, Xinglong Observatory, National Astronomical Observatory of Chinese Academy of Sciences.



#### \*Zyla HF - Fiber Optic coupled Zyla

Check out Andor's new Zyla HF Fiber Optic coupled camera, superb for fast indirect X-Ray applications such as tomography or non-destructive testing.

# ROLLING & GLOBAL SHUTTER

The **Zyla 5.5** uniquely offers both Rolling and *true* Global Shutter exposure modes. This provides superior application and synchronization flexibility and the ability, through global exposure, to closely emulate the familiar ‘Snapshot’ exposure mechanism of interline CCDs.

## Key Benefits of *True* Global Exposure

Global exposure in particular is viewed as an important mode for the biologist, as its benefits are deeply synergistic with the core imaging requirements of live cell microscopy.

- **NO Spatial Distortion** – avoiding the spatial distortion risk of rolling exposure
- Recommended for **3D / 4D microscopy** – Tight syncing to **peripheral switching devices**
- **Higher Signal to Noise** due to **reduced dead time** – the entire exposure cycle can be used
- **Simplicity** – all the benefits of an ‘**interline exposure mode**’
- **Continuous or Pulsed** light sources
- **Sub-microsecond inter-frame gaps** in PIV applications

## ‘Simulated’ Global Exposure in Zyla 4.2 PLUS

[Click here](#) to read more about this mode and other Frequently Asked Questions on Rolling and Global Exposure modes.

## Rolling & Global Shutter Mechanisms


Rolling and *true* Global Shutter modes describe two distinct types of exposure and readout sequence.

In rolling shutter, available in Zyla 4.2 PLUS and Zyla 5.5, different lines of the array are exposed at different times as the read out ‘wave’ sweeps through the sensor. 10 ms is required at the start to ‘activate’ the sensor to expose, and then 10 ms is required at the end to readout the sensor. Use when not synchronizing to peripheral devices AND only when there is a minimal risk of spatial distortion from moving samples.

In *true* global shutter, available in Zyla 5.5, each pixel in the sensor begins the exposure simultaneously and ends the exposure simultaneously. This provides a true ‘Snapshot’ exposure capability for moving samples that is both ‘photon-efficient’ and easy to synchronize to, especially useful for 3D / 4D microscopy. Zyla 4.2 PLUS, while utilizing a rolling shutter sensor, offers a *Simulated* Global Exposure mechanism to overcome risk of spatial distortion. This mechanism is more elaborate and less photon/time efficient than *true* Global Shutter.

[Click here](#) to read more about Rolling and Global shutter modes on our Zyla camera.

### Global Shutter exposure and readout (single scan)

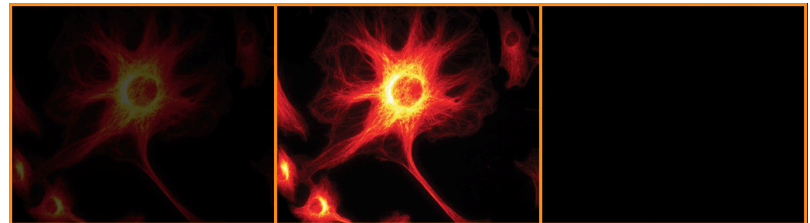


Exposure Start

Exposure

Exposure End

### Rolling Shutter exposure and readout (single scan)

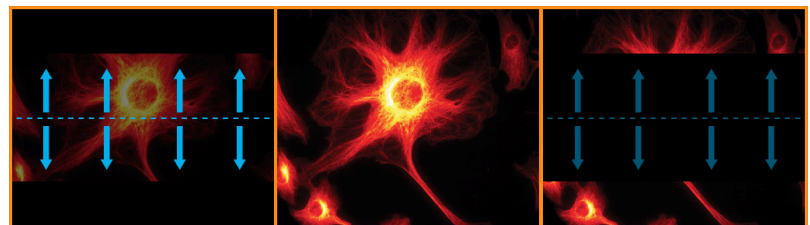

Exposure Start

Exposure

Readout

For further information of Rolling and Global Shutter, please access the following technical notes through the Andor Learning Centre: 1) Rolling and Global Shutter 2) Synchronizing to Rolling and Global Shutter sCMOS cameras

## NEW LightScanPLUS for Zyla 4.2 PLUS

### Key Benefits:

#### FlexiScan


- independent control of pixel row height, scan speed and exposure time
- optimize signal to noise AND confocality concurrently
- multiple read-out options
- scan synchronization output for easy syncing to laser beam

#### CycleMax


- maximum frame rates with reduced dead-time, no need to reset light sheet for each alternate frame

[Click here](#) to find out more about LightScan PLUS.


### Simultaneous Readout Direction Options

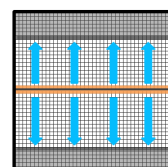


Center Out

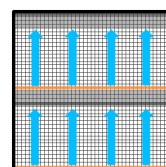

Bottom Up

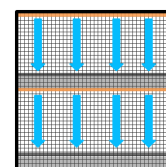


Top Down

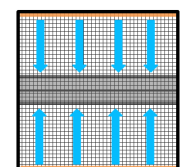


Outside In

### Sequential Readout Direction Options

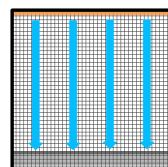

Top Down

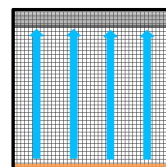

Bottom Up




# sCMOS or EMCCD?

Since the market introduction of sCMOS technology by Andor, the question of the performance comparison against the more established Electron Multiplying CCD (EMCCD) has been common.


Being a very fast, low noise technology, sCMOS does hold some potential to offer an alternative technology to these single photon sensitive detectors across some applications and techniques, including to an extent, super-resolution microscopy and TIRF microscopy. Whilst the read noise of sCMOS is very low compared to CCDs, EMCCD technology holds the distinct advantage of being able to practically eliminate read noise, rendering them single photon sensitive.

After the first few years of sCMOS being in the market, we are concluding that the primary applications for which EMCCDs were originally purchased, such as single molecule detection and low light spinning disk confocal microscopy, are continuing to benefit from this ultrasensitive technology. EMCCDs offer a raw sensitivity that cannot be surpassed in the very low light regime. However, EMCCDs remain relatively expensive, so they will always be considered a more selective, 'high-end' solution.

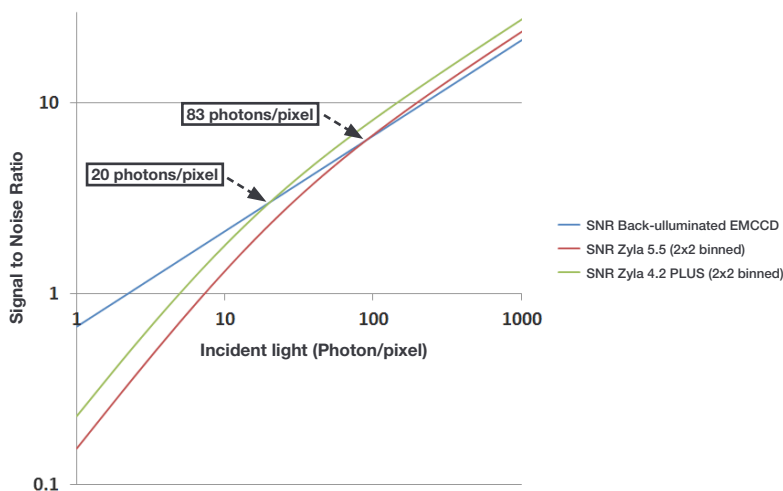

**Figure 1**

Plot of Signal to Noise Ratio versus Incident Photon Intensity, comparing back-illuminated EMCCD iXon 888 (13  $\mu\text{m}$  pixel size) to 2x2 binned Zyla sCMOS cameras (13  $\mu\text{m}$  pixel size after binning). An average QE value for each sensor between 500-750 nm was used.

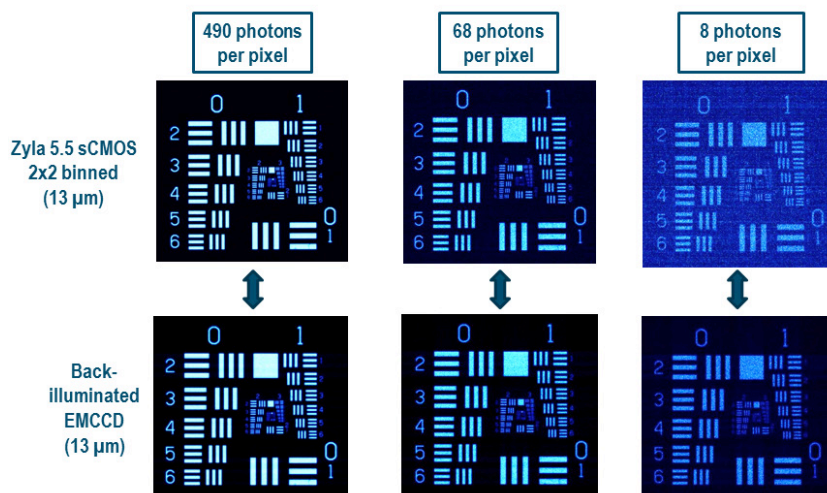

**Figure 2**

Images at a range of incident light intensity, acquired using back-illuminated EMCCD iXon 888 and Zyla 5.5 sCMOS cameras (2x2 binned pixels). At low light intensities, the Signal to Noise Ratio advantage of the EMCCD is apparent.

# TECHNICAL SPECIFICATIONS

## MODEL SPECIFIC SPECIFICATIONS<sup>\*1</sup>

Model	Zyla 5.5	Zyla 4.2 PLUS
<b>Sensor type</b>	Front Illuminated Scientific CMOS	Front Illuminated Scientific CMOS
<b>Active pixels (W x H)</b>	2560 x 2160 (5.5 Megapixel)	2048 x 2048 (4.2 Megapixel)
<b>Sensor size</b>	16.6 x 14.0 mm 21.8 mm diagonal	13.3 x 13.3 mm 18.8 mm diagonal
<b>Pixel readout rate (MHz)</b>	200 (100 MHz x 2 sensor halves) 560 (280 MHz x 2 sensor halves)	Slow Read 216 (108 MHz x 2 sensor halves) Fast Read 540 (270 MHz x 2 sensor halves)
<b>Read noise (e<sup>-</sup>) Median [rms] <sup>*2</sup></b>	<b>Rolling Shutter</b>	<b>Rolling Shutter</b>
	<b>Global Shutter</b>	<b>Global Shutter</b>
	@ 200 MHz 1.2 [1.7] @ 560 MHz 1.45 [1.8]	@ 216 MHz 0.90 [1.4] @ 540 MHz 1.10 [1.6]
<b>Maximum Quantum Efficiency <sup>*3</sup></b>	60%	82%
<b>Sensor Operating Temperature</b>		
<b>Air cooled</b>	0°C (up to 30°C ambient)	0°C (up to 27°C ambient)
<b>Water cooled</b>	-10°C*	-10°C*
<b>Dark current, e<sup>-</sup>/pixel/sec @ min temp <sup>*4</sup></b>		
<b>Air cooled</b>	0.14	0.14
<b>Water cooled</b>	0.04	0.04
<b>Readout modes</b>	Rolling Shutter and True Global Shutter (Snapshot)	Rolling Shutter and Global Clear <sup>*8</sup>
<b>Maximum dynamic range</b>	25,000:1	33,000:1
<b>Photon Response Non-Uniformity (PRNU)</b>	< 0.5%	< 0.1%
<b>Pre-defined Region of Interest (ROI)</b>	2048 x 2048, 1920 x 1080, 1392 x 1040, 512 x 512, 128 x 128	1920 x 1080, 1392 x 1040, 512 x 512, 128 x 128
<b>User defined ROI granularity</b>	1 pixel <sup>**</sup>	
<b>Data range</b>	12-bit (fastest USB 3.0 speeds) and 16-bit (maximum dynamic range)	
<b>Interface options</b>	USB 3.0 <sup>*9</sup> Camera Link 10-tap	

\* Cooling temperature must be above the dew point

\*\* Minimum ROI size possible: 16 x 12 in 12-bit mode and 12 x 12 in 16-bit mode

## GENERAL SPECIFICATIONS<sup>\*1</sup>


<b>Pixel size (W x H)</b>	6.5 µm
<b>Pixel well depth (e<sup>-</sup>)</b>	30,000
<b>Linearity (% , maximum) <sup>*5</sup></b>	Better than 99.8%
<b>MTF (Nyquist @ 555 nm)</b>	45%
<b>Pixel binning</b>	Hardware binning: 2 x 2, 3 x 3, 4 x 4, 8 x 8
<b>Anti-blooming factor</b>	x 10,000
<b>I/O</b>	External Trigger, Fire, Fire n, Fire All, Fire Any, Arm
<b>Trigger Modes</b>	Internal, External, External Start, External Exposure, Software Trigger
<b>Software Exposure Events<sup>*6</sup></b>	Start exposure - End exposure (row 1), Start exposure - End exposure (row n)
<b>Hardware timestamp accuracy</b>	25 ns
<b>Internal memory</b>	1 GB




## FRAME RATE TABLE - 12-BIT (16-BIT)<sup>\*7</sup>

Array Size	Zyla 5.5 USB 3.0		Zyla 5.5 10-tap		Zyla 4.2 PLUS 10-tap	Zyla 4.2 PLUS USB 3.0
	Rolling Shutter	Global Shutter	Rolling Shutter	Global Shutter	Rolling Shutter	Rolling Shutter
2560 x 2160	40 (30)	40 (30)	100 (75)	49 (49)	-	-
2048 x 2048	53 (40)	52 (39)	105 (98)	52 (52)	101 (101)	53 (40)
1920 x 1080	107 (80)	98 (80)	200 (200)	97 (97)	192 (192)	107 (80)
512 x 512	422 (422)	201 (201)	422 (422)	201 (201)	406 (406)	406 (406)
128 x 128	1691 (1691)	716 (716)	1691 (1691)	716 (716)	1627 (1627)	1627 (1627)
2048 x 8 (FCS mode)	27057 (27057)	4008 (4008)	27057 (27057)	4008 (4008)	26041 (26041)	26041 (26041)

## QUANTUM EFFICIENCY (QE) CURVE<sup>\*3</sup>

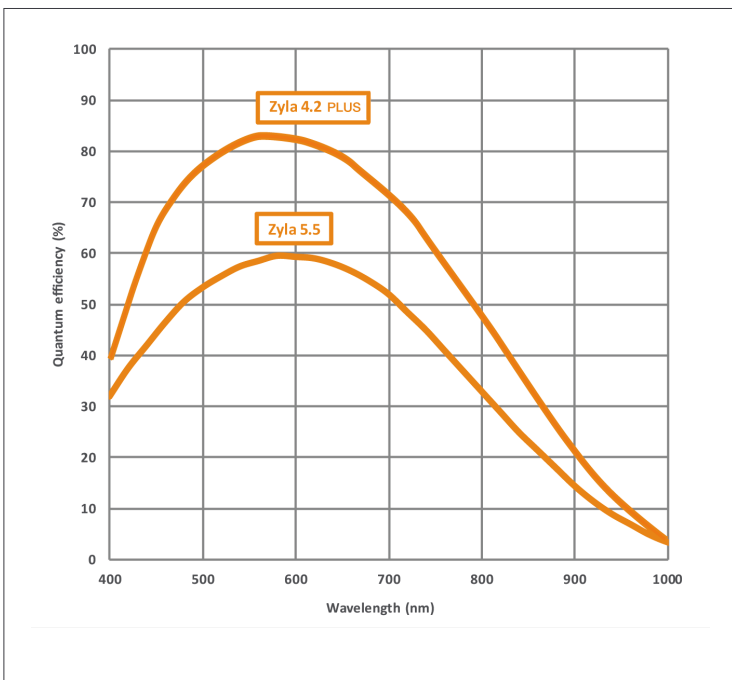

## QE VS. FLUOROPHORE EMISSIONS

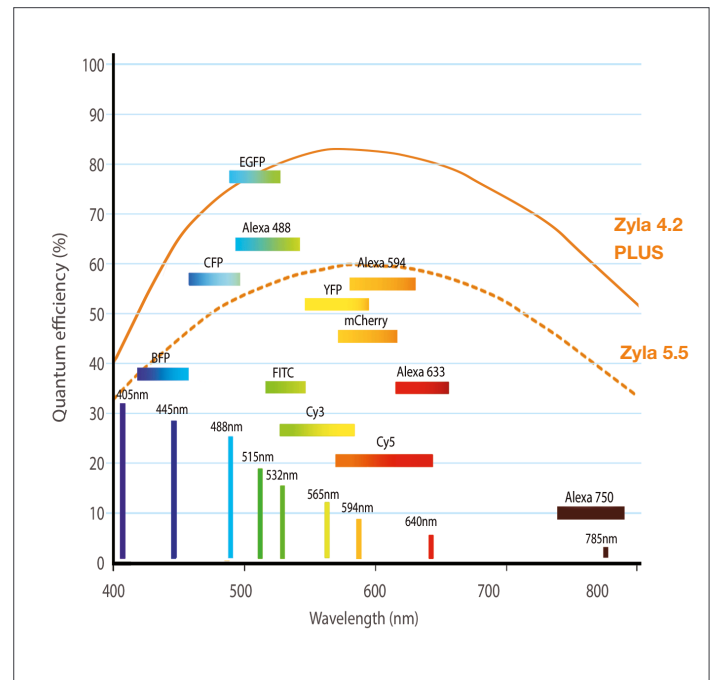

## GPU Express




The Andor GPU Express library has been created to simplify and optimize data transfers from camera to a CUDA-enabled NVidia Graphical Processing Unit (GPU) card to facilitate accelerated GPU processing as part of the acquisition pipeline. GPU Express integrates easily with SDK3 for Windows, providing a user-friendly but powerful solution for management of high bandwidth data flow challenges; ideal for data intensive applications such as Light Sheet Microscopy, Super-Resolution Microscopy and Adaptive Optics.

- Enhanced convenience, afforded by simple, optimized GPU data management
- Optimal data throughput
- Superb, easily accessible documentation and examples.

## CREATING THE OPTIMUM PRODUCT FOR YOU

### Step 1. Select the camera type

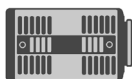


Camera Type

Description	Code
ZYLA 4.2 PLUS, 4.2 Megapixel, Rolling shutter, 100 fps, Camera Link 10-tap	<b>ZYLA-4.2P-CL10</b>
ZYLA 4.2 PLUS, 4.2 Megapixel, Rolling shutter, 53 fps, USB 3.0	<b>ZYLA-4.2P-USB3</b>
ZYLA 5.5, 5.5 Megapixel, Rolling and Global shutter, 100 fps, Camera Link 10-tap	<b>ZYLA-5.5-CL10</b>
ZYLA 5.5, 5.5 Megapixel, Rolling and Global shutter, 40 fps, USB 3.0	<b>ZYLA-5.5-USB3</b>

For water cooled option, add **-W** to your selected camera code

### Step 2. Select the required accessories

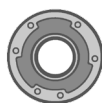

Accessories

Description	Order Code
CS-mount adapter	<b>ACC-MEC-05609</b>
F-mount adapter	<b>ACM-05574</b>
Auto extension tubes (set of 3) for C-mount	<b>OA-ECMT</b>
Auto extension tubes (set of 3) for Nikon F	<b>OA-ENAF</b>
Re-circulator for enhanced cooling performance	<b>XW-RECR</b>
Oasis 160 Ultra compact chiller unit	<b>ACC-XW-CHIL-160</b>
5 meter Camera Link connector cable. Order x2 if using with Zyla Camera Link 10-tap models.	<b>ACC-ASE-02992</b>
10 meter active Camera Link connector cable, including power supply. For use with Zyla 10-tap Camera Link models.	<b>ACC-ASE-06962</b>
30 meter fibre-optic extender solution for use with Zyla Camera Link 10-tap models.	<b>ACC-ZYLFOX-10TAP-30M</b>
100 meter fibre-optic extender solution for use with Zyla Camera Link 10-tap models.	<b>ACC-ZYLFOX-10TAP-100</b>
15 meter active USB 3.0 connector cable (power supply not required). For use with Zyla USB 3.0 models.	<b>ACC-ASE-06887</b>
50m fibre optic USB 3.0 extender solution including power supply. For use with Zyla USB 3.0 models.	<b>ACC-ASE-08762</b>
100m fibre optic USB 3.0 extender solution including power supply. For use with Zyla USB 3.0 models.	<b>ACC-ASE-07860</b>
PC Workstation for up to 100 fps continuous spooling to hard drives, acquiring up to 120,000 12-bit full resolution images: Dell T7910XL, 2.6 GHz Eight Core, 8 GB RAM, 4 x 250GB SSD hard drive configured in RAID 0.	<b>WKST-1 WIN</b>
PC Workstation for up to 100 fps continuous spooling to RAM, acquiring up to 6,000 12-bit full resolution images: Dell T5810, 3.5 GHz Quad Core, 64 GB RAM.	<b>WKST-3 WIN</b>

For further information on PC workstations for Zyla, please refer to the technical note [PC Specifications for sCMOS](#)

### Step 3. Select the required software



Software

**The Zyla also requires at least one of the following software options:**

**Solis Imaging** A 32-bit and fully 64-bit enabled application for Windows (7 and 8) offering rich functionality for data acquisition and processing. AndorBasic provides macro language control of data acquisition, processing, display and export.

**Andor iQ** A comprehensive multi-dimensional imaging software package. Offers tight synchronization of the camera with a comprehensive range of microscopy hardware, along with comprehensive rendering and analysis functionality. Modular architecture for best price/performance package on the market.

**Andor SDK3** A software development kit that allows you to control the Andor sCMOS cameras from your own application. Available as 32 and 64-bit libraries for Windows (7 and 8) and Linux. Compatible with C/C++, LabView and Matlab.




**GPU Express** Andor GPU Express library has been created to simplify and optimize data transfers from camera to a CUDA-enabled NVidia Graphical Processing Unit (GPU) card to facilitate accelerated GPU processing as part of the acquisition pipeline. Integrates easily with Andor SDK3 for Windows.

**Third party software compatibility**

Drivers are available so that the Zyla can be operated through a large variety of third party imaging packages. See Andor web site for detail: [andor.com/software/](http://andor.com/software/)

## PRODUCT DRAWINGS


Dimensions in mm [inches]

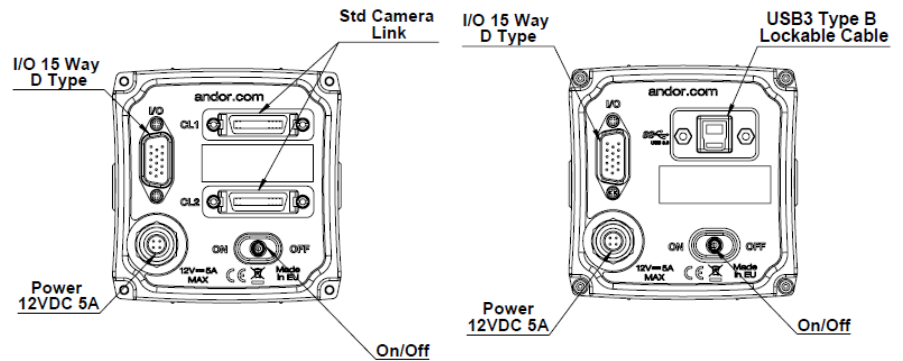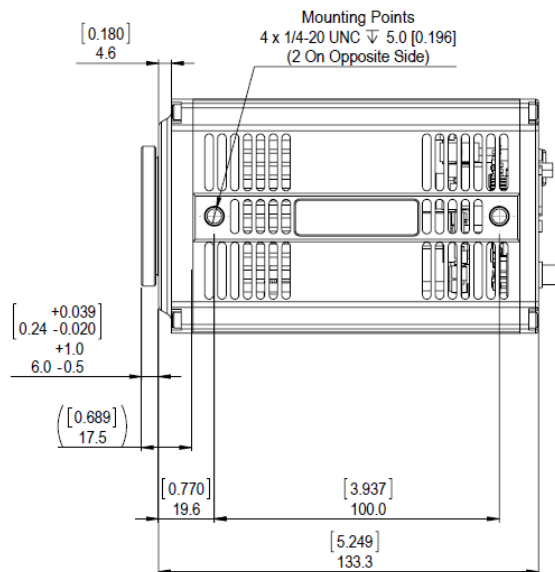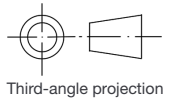

**Note:** Please leave 70mm minimum for the bend radius of the USB 3.0 cable.

**Weight:** 1,000 g [2 lbs 3 oz]

Product drawings of the water cooled Zyla can be found at <http://www.andor.com/watercooledzyla>

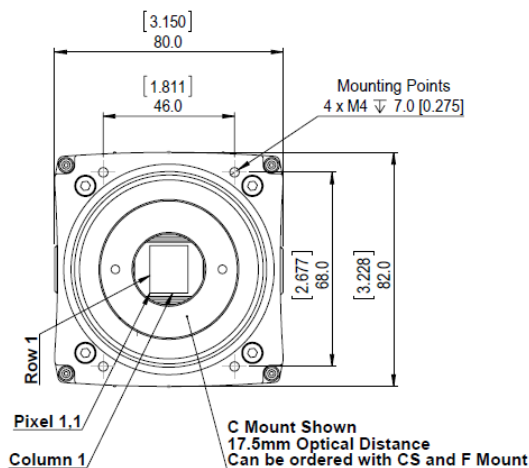

## CONNECTING TO THE ZYLA

### Camera Control

Connector type: 3 meter Camera Link 10-tap connectors or USB 3.0. (Longer lengths available as accessories).

### TTL / Logic

Connector type: 15 way D Type with TTL I/Os for External Trigger, Frame Readout and Fire Pulse

### REGULATORY COMPLIANCE

- RoHS compliant
- EU EMC Directive
- EU LV Directive
- IEC 61010-1 CB Scheme

### EXTERNAL POWER SUPPLY COMPLIANCE

- UL-certified for Canada and USA
- Japanese PSE Mark

### POWER SUPPLY REQUIREMENTS

- Power: +12 VDC  $\pm$  5% @ 5A
- Ripple: 200 mV peak-peak 0 - 20 MHz
- 100 - 240 VAC 50/60 Hz external power supply
- Power Consumption: 12V @ 5A Max, 12V @ 2.5A Nominal

## 15-WAY D-TYPE PINOUTS

1	ARM	Output
2	Aux_Out_1*	Output
3	FIRE row n	Output
4	FIRE row 1	Output
5	Aux_Out_2	Output
6	Ground	GND
7	External Trigger	Input
8	Spare Input	Input
9	Reserved	N/A
10	Reserved	N/A
11	Reserved	N/A
12	Reserved	N/A
13	Reserved	N/A
14	Reserved	N/A
15	Reserved	N/A

\* Aux\_Out\_1 is configurable as Fire, Fire n, Fire All or Fire Any. Refer to the Zyla hardware manual.



# ORDER TODAY

Need more information? At Andor we are committed to finding the correct solution for you. With a dedicated team of technical advisors, we are able to offer you one-to-one guidance and technical support on all Andor products. For a full listing of our regional sales offices, please see: [andor.com/contact](http://andor.com/contact)

Our regional headquarters are:

### EUROPE

Belfast, Northern Ireland  
Phone +44 (28) 9023 7126  
Fax +44 (28) 9031 0792

### JAPAN

Tokyo  
Phone +81 (3) 6732 8968  
Fax +81 (3) 6732 8939

### NORTH AMERICA

Connecticut, USA  
Phone +1 (860) 290 9211  
Fax +1 (860) 290 9566

### CHINA

Beijing  
Phone +86 (10) 8271 9066  
Fax +86 (10) 8271 9055

### ITEMS SHIPPED WITH YOUR CAMERA

**For Camera Link 10-Tap Models:** 1 x Camera Link Card and 2 x 3 meter connector cables.

**For USB 3.0 models:** 1 x USB 3.0 PCIe Card and 1 x 3 meter USB 3.0 cable (Type A to B)  
1 x Power supply with mains cable  
1 x 7-way Multi I/O timing cable, offering Fire, External Trigger and Arm (3 meter)  
1 x Quick Start Guide  
1 x CD containing Andor user guides

### FOOTNOTES:

- Specifications are subject to change without notice
- Figures are typical unless otherwise stated.
  - Readout noise is for the entire system and is taken as a median over the sensor area excluding any regions of blemishes. It is a combination of sensor readout noise and A/D noise.
  - Quantum efficiency of the sensor at 20°C as supplied by the manufacturer.
  - Dark current measurement is taken as a median over the sensor area excluding any regions of blemishes.
  - Linearity is measured from a plot of Signal vs. Exposure Time over the full dynamic range.
  - Software Exposure Events provide rapid software notification (SDK only) of the start and end of acquisition, useful for tight synchronization to moving peripheral devices e.g. Z-stage.
  - The maximum frames/s table for Zyla indicate the maximum speed at which the device can acquire images in a standard system at full frame and also a range of sub-array size, for both rolling and global shutter read modes (Zyla 5.5), 12-bit single amplifier (rates also apply to dual amplifier 16-bit for Zyla 4.2). Note that the write speed of the PC hard drive can impose a further restriction to achieving sustained kinetic series acquisition.
  - 'Global Clear' is an optional keep clean mechanism that can be implemented in rolling shutter mode, which purges charge from all rows of the sensor simultaneously, at the exposure start. The exposure end is still rolling shutter. It can be used alongside the Fire All output of the camera and a pulsed light source to simulate Global Exposure mechanism, albeit less efficiently than the true Global Shutter exposure mode of Zyla 5.5. Furthermore Global Clear differs from true Global Shutter in that it can only be used in 'non-overlap' readout mode, i.e. sequential exposure and readout phases rather than simultaneous.
  - Zyla USB 3.0 models should work with any modern USB 3.0 enabled PC/laptop (provided hard drives or RAM is sufficient to support data rates) as every USB 3.0 port should have its own host controller. Zyla USB 3.0 models also ship with a USB 3.0 PCI card as a means to add a USB 3.0 port to an older PC, or as a diagnostic aid to interoperability issues or to ensure maximum speed.



### MINIMUM COMPUTER REQUIREMENTS:

- 2.68 GHz Quad Core
- 4GB RAM (increase RAM if to be used for continuous data spooling)
- Hard Drive:  
Minimum 450 MB/s continuous write for USB 3.0 models  
Minimum 850 MB/s continuous write for Camera Link 10-tap models
- PCI Express x4 only for USB 3.0 models
- PCI Express x8 or greater for Camera Link 10-tap models
- Windows (7 or 8) or Linux
- \* See technical note entitled: 'PC Specifications for sCMOS'
- \*\* Note, Andor supply PC workstations for Zyla, see page 10.

### Operating and Storage Conditions

- Operating Temperature:  
Zyla 5.5: 0°C to 30°C ambient  
Zyla 4.2: 0°C to 27°C ambient
- Relative Humidity: < 70% (non-condensing)
- Storage Temperature: -10°C to 50°C

### Power Requirements

- Please refer to page 11



Windows is a registered trademark of Microsoft Corporation